



Key International Inc. 480 Rt. 9 Englishtown, NJ 07726  
(732) 536-9700 Phone (732) 972-2630 Fax [www.keyinternational.com](http://www.keyinternational.com)

### **HT-300/500 and 300S/500S Quick Start Guide**

Listed below please find the basic operational procedure for the Key Tablet Hardness Tester:

- 1.) Apply power and turn unit on via Power Switch. Printer will print a row of zeroes.
- 2.) Press the X,O button to turn on the Statistics function. LED will illuminate.
- 3.) Set the Continuous restart value using the CONT button, if desired.
- 4.) Place the first sample on the pedestal, then close isolation cover.
- 5.) Press START. Motor will initiate test cycle.
- 6.) When the sample breaks, a result will be printed. Brush off tablet pieces.
- 7.) Place next sample onto pedestal and wait for Automatic Restart or press START.
- 8.) Continue following procedure from step 6, until all samples have been tested.
- 9.) Press STOP to end testing.
- 10.) Press X,O to print out statistical report.

END