



**Key International Inc. —**  
**4 Corporate Dr.**  
**Cranbury, NJ 08512**  
**Phone: 609-619-3685 Fax: 609-619-3686**  
**www.keyinternational.com**  
**info@keyinternational.com**

---

## **HT-300/500 - Unit Selection Procedure**

**Valid for units with Serial Numbers 4500 and above**

- 1.) Press and hold STOP, then press START. Release both keys. A message will print - "PTB 301 xx.xx/x".
- 2.) Press and hold START while watching the display. It will scroll through a series of numbers – 0 1 2 3.
- 3.) Release START when the desired unit is reached as shown below:

0	=	Newtons	(N)	
1	=	Kiloponds	(KP)	
2	=	Strong Cobbs	(SC)	
3	=	Digits	(D)	(Used only for factory testing)

- 4.) Once a unit is selected, it will be printed on the printer.
- 5.) Press STOP. A confirming row of zeroes will print.